


# pCLIP<sub>f</sub>-H2B Control Plasmid



N9218S 006120615061

## N9218S

**20 µg** Lot: 0061206  
**Store at: -20°C** Exp: 6/15

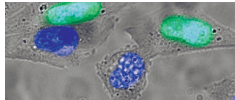

Live CHO-K1 cells transiently transfected with pCLIP<sub>f</sub>-H2B. Cells were labeled with CLIP-Cell™ 505 (green) for 30 minutes and counterstained with Hoechst 33342 (blue).

### Introduction

This control plasmid contains the gene encoding the Histone H2B protein cloned upstream of the CLIP<sub>f</sub> coding sequence in pCLIP<sub>f</sub> as a fusion to the N-terminus of the CLIP-tag. Histone H2B is a member of the core histones that package DNA in the nucleus. The H2B-CLIP<sub>f</sub> fusion protein gives nuclear fluorescence when labeled with CLIP-Cell™ substrates. The full sequence and map for pCLIP<sub>f</sub>-H2B can be downloaded at [www.neb.com](http://www.neb.com).

The CLIP-tag is a novel tool for protein research, allowing the specific, covalent attachment of virtually any molecule to a protein of interest. The CLIP-tag is a small polypeptide based on human O<sup>6</sup>-alkylguanine-DNA-alkyltransferase (AGT). CLIP-tag substrates are derivatives of benzylcytosine (BC). In the labeling reaction, the substituted benzyl group of the substrate is covalently attached to the reactive cysteine of CLIP-tag forming a stable thioether link. Although CLIP-tag is based on the same protein as SNAP-tag®, the benzylcytosine substrates form a separate class of substrates, different from the benzylguanine substrates recognized by SNAP-tag. CLIP-tag and SNAP-tag can be used for orthogonal simultaneous labeling.

pCLIP<sub>f</sub> contains an improved version of CLIP-tag, termed CLIP<sub>f</sub>. CLIP<sub>f</sub> displays faster kinetics in *in vitro* labeling and fast, specific and efficient labeling in live and fixed cell applications, thereby rendering it a desired research tool for analysis of protein dynamics.

There are two steps to using this system: sub-cloning and expression of the protein of interest as a CLIP<sub>f</sub> fusion, and labeling of the fusion with the CLIP-tag substrate of choice. Expression of the CLIP<sub>f</sub>-H2B fusion protein is described in this document. The labeling of fusion proteins with CLIP-tag substrates is described in the instructions supplied with CLIP-tag substrates.

### Materials Required but not Supplied:

Tissue culture media and reagents  
Mammalian cell lines  
Transfection reagents  
CLIP-tag substrates

### Storage

pCLIP<sub>f</sub>-H2B is supplied in TE buffer (10 mM Tris-HCl, pH 8.0, 1 mM EDTA) at a concentration of 0.5 µg/µl. Plasmid solutions can be stored at 4°C for up to one week. For long-term storage -20°C is recommended.

### Expression of CLIP<sub>f</sub> Fusions


#### Transient Expression

Expression of the fusion protein cloned in pCLIP<sub>f</sub>-H2B can be achieved by transiently transfecting cells in culture with standard transfection protocols. The appropriate reagent and time to permit adequate expression must be empirically determined. pCLIP<sub>f</sub>-H2B has performed well in stable and transient transfection of CHO-K1, COS-7, U-2 OS and NIH 3T3 cells. Note that the intensity of the fluorescence may vary depending on cell line and labeling substrate used.

We recommend using TransPass D2 (NEB #M2554) in combination with TransPass V (NEB #M2561) or Roche's FuGENE® 6 Transfection Reagent for both transient and stable transfections.

#### Stable Expression

pCLIP<sub>f</sub>-H2B can be transfected as described above for transient transfection or by other standard transfection methods. Twenty four to 48 hours after transfection, begin selecting mammalian cultures in 600–1,200 µg/ml G418 (geneticin) depending on the cell line. It is recommended that a kill curve be established for each cell line to determine optimal selection conditions. After 8–12 days of continuous selection, stable colonies will become visible. It is possible to use pools of stable cell populations for initial cell labeling to test for the presence of CLIP-tag expression. In addition, clonal cell lines can be isolated and characterized if desired.

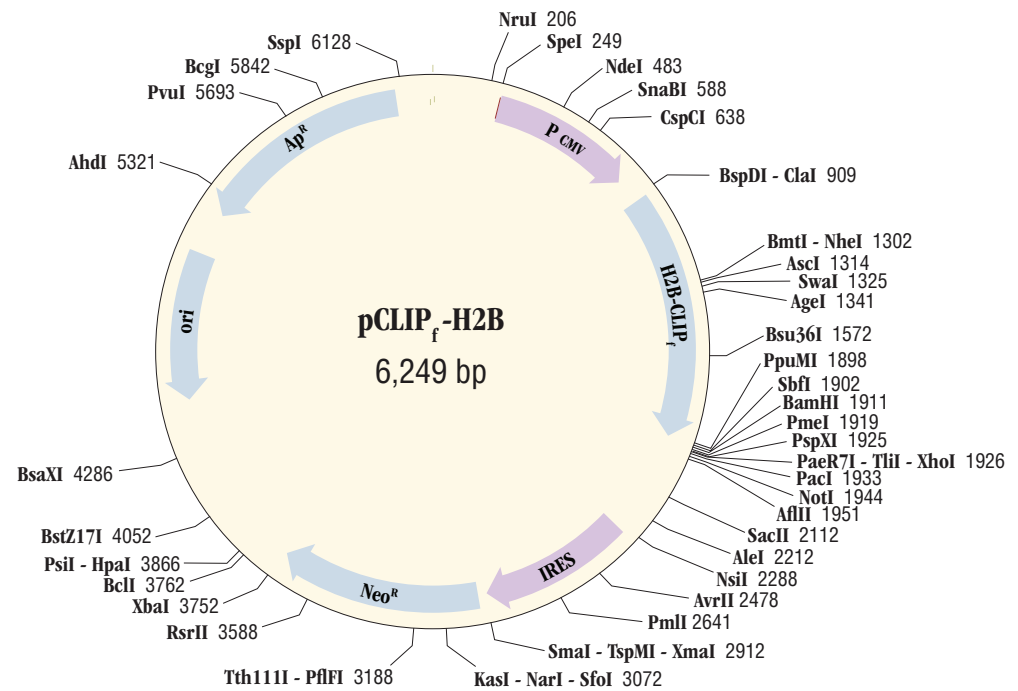

### Troubleshooting

#### Expression

In general, we have not experienced problems expressing H2B-CLIP<sub>f</sub> from the pCLIP<sub>f</sub>-H2B plasmid. Labeling of transfected cells with a fluorescent CLIP-Cell substrate should show strong nuclear fluorescence. In most instances, difficulties in expression can be resolved by altering the transfection protocol.

The CMV promoter is covered under U.S. Patent No. 5,385,839 and its use is permitted for research purposes only. Any other use of the CMV promoter requires a license from the University of Iowa Research Foundation, 214 Technology Innovation Center, Iowa City, IA 52242.

**Notice to Buyer/User:** The Buyer/User has a non-exclusive license to use this system or any component thereof for **RESEARCH AND DEVELOPMENT PURPOSES ONLY**. Commercial use of this system or any components thereof requires a license from New England Biolabs, Inc., 240 County Road Ipswich, MA 01938. For detailed information, see: [www.neb.com/cia/legal](http://www.neb.com/cia/legal).

**For Research Use Only:** No other use is authorized, including without limitation Commercial Use. Commercial Use means any and all uses of the product and derivatives by a party for monetary or other consideration including, but not limited to: (1) product manufacture; (2) providing of a service, information or data; (3) resale of the product, derivatives or components; (4) use of the product, derivatives or components by therapeutic, diagnostic or prophylactic clinical or commercial purposes. Information regarding any other license may be obtained by contacting Business Development and Licensing, New England Biolabs, Inc., 240 County Road, Ipswich, MA 01938. Tel: 978-927-5054, Fax: 978-921-1350.

FuGENE® is a registered trademark of Roche.